

## Running a Spectrum Unlocked, with a Z-Gradient

1. Lock, shim, and run a standard spectrum, as usual.
2. Turn the lock and lock-sweep off:  
In the *Lock/Level* tab of the `BSMS` control panel, unclick the `lock` and `sweep` buttons.
3. If running with a gradient,
  - (a) change `pulprog` to `zg30Grad`
  - (b) set `gpz0` to the desired z-gradient strength. (0 – 100% of maximum; 0.1 – 1 work well)
4. Run spectrum (`zg`) as usual.
5. When done, don't forget
  - to turn the lock-sweep and lock back on, in the `BSMS` panel.
  - to reset `gpz0` to 0.